



Automated Sample Preparation

# **DYNAMIC HEADSPACE SYSTEM +**


**DHS+**

Fully Automated VOC Extraction

# Dynamic enrichment for maximum sensitivity in GC-MS analysis

GERSTEL Dynamic Headspace<sup>+</sup> (DHS<sup>+</sup>) is a fully automated, dynamic headspace technique for highly efficient enrichment of volatile and semi-volatile compounds from complex sample matrices – even at low extraction temperatures.

The sample is tempered, agitated, or cooled in a sealed vial. Volatile compounds are continuously purged from the headspace, enriched on an adsorbent tube, and then automatically thermally desorbed, refocused, and transferred Quantitatively onto the GC column.

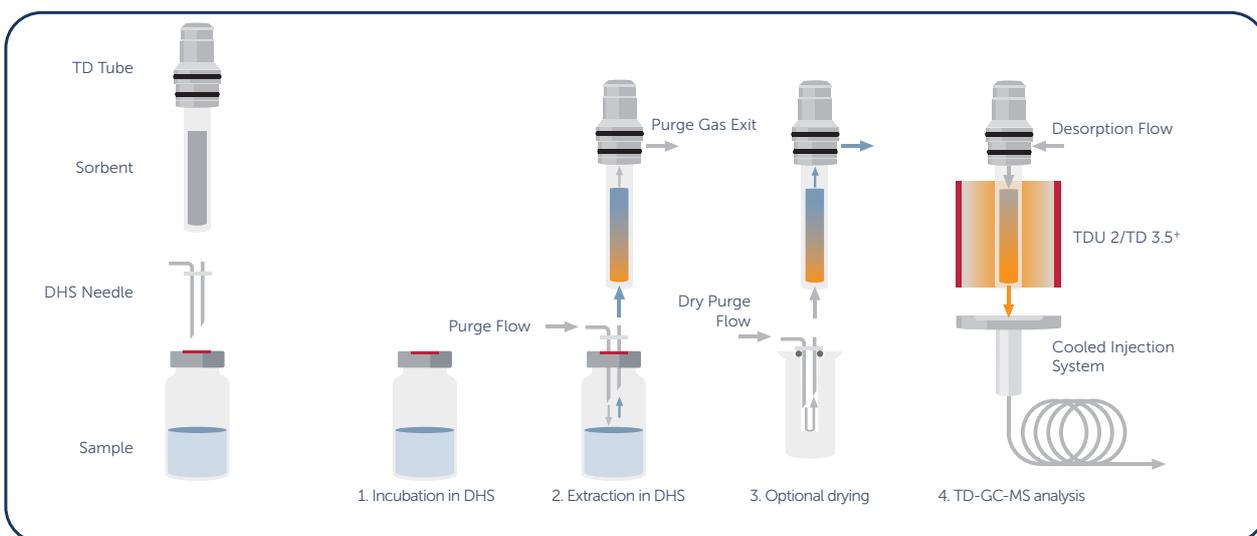

**Schematic representation of the DHS<sup>+</sup> workflow:** Continuous extraction of the analytes from the headspace, enrichment on the adsorbent tube, thermal desorption and refocusing prior to GC-MS analysis.


## What makes DHS<sup>+</sup> special

- High sensitivity even at low extraction temperatures
- Enrichment until complete sample depletion
- Very low detection limits and large linear range
- Wide extraction bandwidth: VVOC to SVOC and polar to nonpolar
- Independent of equilibrium settings, as with SPME

## Was DHS<sup>+</sup> besonders macht


- Continuous extraction with controlled inert gas flow
- Multi-sorbent trapping for maximum analyte coverage
- Fractionated DHS. multiple extractions from one sample
- Desorption of multiple traps / extracts into one GC run
- Optional: Cryogen-free focusing via dynamic focusing
- Automatically exchangeable traps to prevent carryover
- Automated monitoring of flow stability and system tightness
- Ensured reproducible recoveries through controlled extraction integrity

**Result:** reproducible quantification even for challenging matrices – from foods and fragrances to environmental and polymer samples.

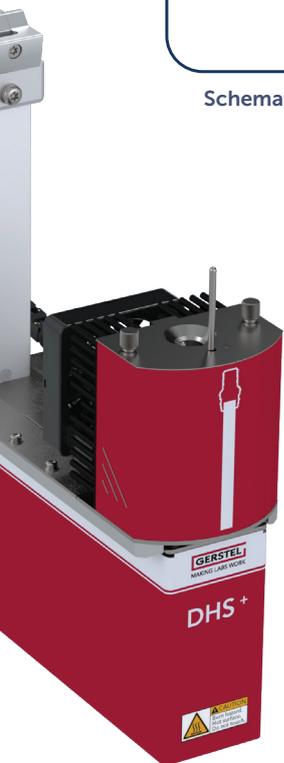


# Matrix-independent VOC analysis with FET and MVM

In the **Full Evaporation Technique (FET)**, the parameters are optimized to completely evaporate analytes using rather gentle temperatures and a huge gas volume in order to resolve matrix effects and then dynamically enriched – regardless of matrix effects or distribution coefficients.

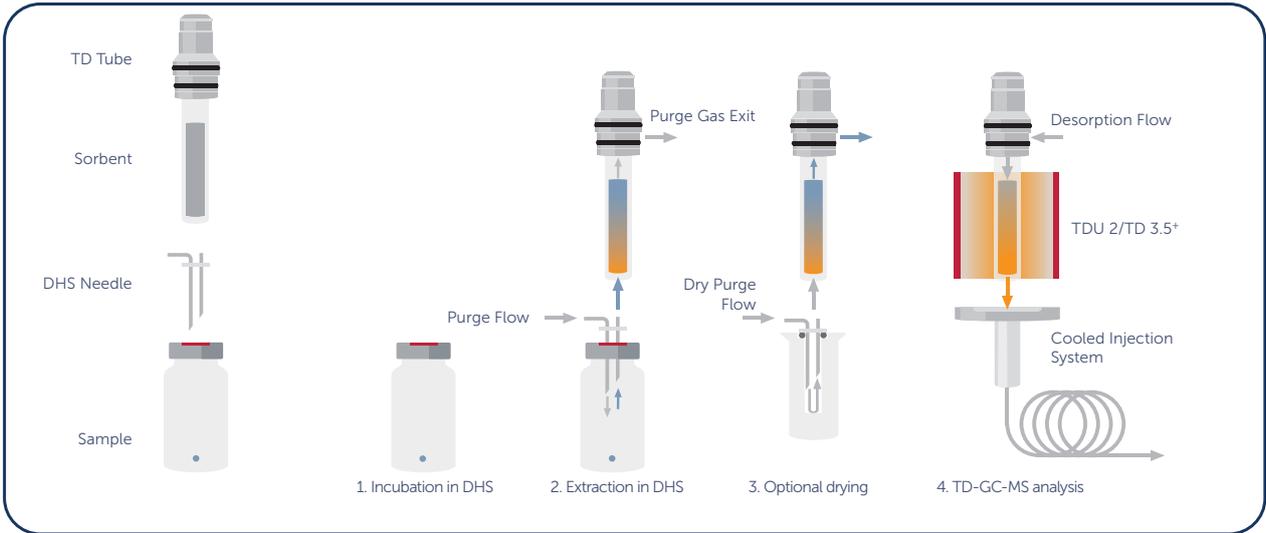

Schematic representation of the Full Evaporation Technique (FET)

The **Multi Volatiles Method (MVM)** is based on the concept of fractionated DHS, where multiple dynamic headspace extracts can be combined to detect compounds with varying volatility. By combining optimized extraction conditions with multi-sorbent trapping, MVM enables the reproducible and matrix-independent analysis of a broad volatility spectrum within a chromatography run. It is ideal for non-target analyses like Beverages or cosmetic products.

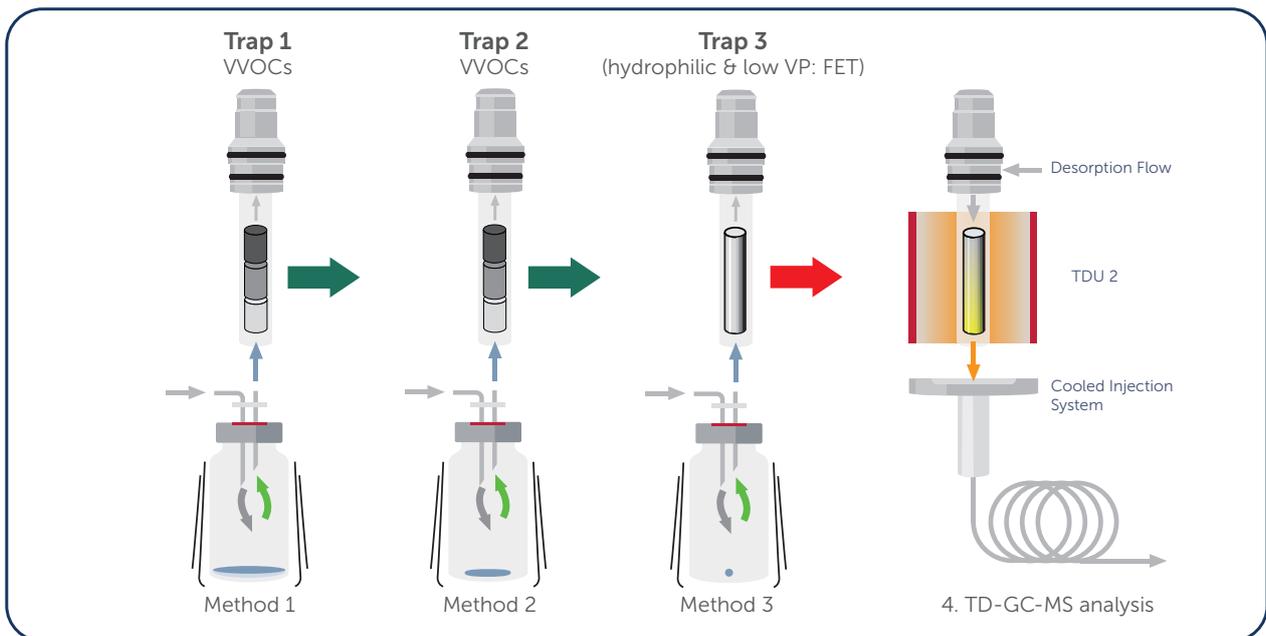

Schematic representation of the Multi Volatile Method (MVM) workflow

# Applications & Benefits

## Typical applications

- Aroma and flavor profiling
- Off-flavor and odor analysis
- Fragrance and market sample analysis
- Extractables & Leachables (E&L)
- Packaging analysis
- Emission analysis of polymers and textile fibers
- Environmental VOC analysis
- Target and non-target screening



## Your benefits in everyday laboratory work

- Ultra-trace VOC analysis with high sensitivity
- Matrix-independent quantification using FET (Full Evaporation Technique) and MVM (Multi Volatiles Method)
- Fully automated workflows – ideal for routine analysis and screening
- Solvent-free, sustainable microextraction technique
- High sample throughput with minimal operator effort
- Flexible quantification strategies via integrated sample preparation e.g. Standard addition

**DHS+** extends the proven DHS principle with additional automation, monitoring, and control functions for increased process reliability, reproducible recoveries, and robust quantification – particularly for demanding and regulated applications.



**GERSTEL GmbH**  
Headquarters

📍 **Eberhard Gerstel Platz 1**  
45473 Mülheim an der Ruhr  
Germany

☎ +49 208 76503-0  
✉ [info@gerstel.com](mailto:info@gerstel.com)  
🌐 [www.gerstel.com](http://www.gerstel.com)

Velaris reserves the right to change the specifications and the appearance of the equipment without further notification.

Version 1.0